

# Add specimen requirement groups0

You may add any number of specimen requirements for a collection event and group them together into groups. To add a specimen requirement group, search for the desired study, using [Search](#) [ #id.elwy3aftu9nv]a [ #id.elwy3aftu9nv]study and navigate to a collection event whose specimen requirements should be grouped.

Study details

Study details Associated study(s)

Alliance study number \* : 10600

Title \* : 10600 - Treatment study

Start date \* : 07/01/2012 14

Author : Admin BMS

Clinical diagnosis : Start typing here

Abnormality of neurotensine secretion (disord  
[M]Dermatofibroma NOS

Submit Add epoch

Save study Synchronize study

Select the collection event from the tree structure

Study details

ADD COLLECTION EVENT

Event label \* : Screening Co-incident event

Clinical status \* : Not Specified

Calendar event point \* : 2.0 Days

Clinical diagnosis \* : [M]Medulloblastoma NOS

SPECIMEN REQUIREMENT GROUPS

Add Delete

Submit Cancel Add Specimen Requirement

Save study Synchronize study

In add collection event page, from the specimen requirement groups section, select *Add* button.

**Study details**

- 10600
  - Pre-Registration
    - Pre-Registration
      - Screening**
        - 5 ml Bone marrow Aspirate (Green Top)
          - Bone Marrow Biopsy Section
        - 2 Bone Marrow Biopsy Sections (for immuno)
        - 6ml Whole Blood (Green top)
        - 6ml Whole Blood (Green top)
        - 6 Unstained Bone Marrow Smears (for cyto)
    - Remission Induction Therapy
    - Midostaurin

### ADD COLLECTION EVENT

Event label \* :  [Co-incident event](#)

Clinical status \* :

Calendar event point \* :  Days

Clinical diagnosis \* :

---

### SPECIMEN REQUIREMENT GROUPS

Group label *	Specimen requirement(s) *
<input type="checkbox"/> 12 ml Whole Blood ( 2 x 6ml Gr	<div style="display: flex; align-items: flex-start;"> <div style="flex: 1;"> <input type="text" value="Please select"/> <div style="border: 1px solid #ccc; padding: 2px; margin-top: 2px;">                     Please select                      5 ml Bone marrow Aspirate (Green Top)                      2 Bone Marrow Biopsy Sections (for immuno)                      6ml Whole Blood (Green top)                      Bone Marrow Biopsy Section                      6ml Whole Blood (Green top)                      6 Unstained Bone Marrow Smears (for cyto)                 </div> </div> <div style="flex: 0 0 60px; text-align: center; margin-left: 5px;"> <input type="button" value="Add &gt;&gt;"/>  <input type="button" value="&lt;&lt; Remove"/> </div> </div>

Enter a Group label and from the dropdown, select the list of specimen requirements that should be grouped together. In the Specimen requirements dropdown, system displays the list of all specimen requirements for the selected collection event. Select the desired specimen requirement and select *Add* button. Now, the added specimen requirement is displayed in selected specimen requirement list. To add another specimen requirement into the group, select it again and select *Add* following which it is added to list of selected ones.

**Study details**

- 10600
  - Pre-Registration
    - Pre-Registration
      - Screening**
        - 5 ml Bone marrow Aspirate (Green Top)
          - Bone Marrow Biopsy Section
        - 2 Bone Marrow Biopsy Sections (for immuno)
        - 6ml Whole Blood (Green top)
        - 6ml Whole Blood (Green top)
        - 6 Unstained Bone Marrow Smears (for cyto)
    - Remission Induction Therapy
    - Midostaurin

### ADD COLLECTION EVENT

Event label \* :  [Co-incident event](#)

Clinical status \* :

Calendar event point \* :  Days

Clinical diagnosis \* :

---

### SPECIMEN REQUIREMENT GROUPS

Group label *	Specimen requirement(s) *
<input type="checkbox"/> 12 ml Whole Blood ( 2 x 6ml Gr	<div style="display: flex; align-items: flex-start;"> <div style="flex: 1;"> <input type="text" value="Please select"/> <div style="border: 1px solid #ccc; padding: 2px; margin-top: 2px;">                     6ml Whole Blood (Green top)                      6ml Whole Blood (Green top)                 </div> </div> <div style="flex: 0 0 60px; text-align: center; margin-left: 5px;"> <input type="button" value="Add &gt;&gt;"/>  <input type="button" value="&lt;&lt; Remove"/> </div> </div>

After adding the desired specimen requirements into the group, select *Submit*. System groups the selected specimen requirements and saves the collection event. You may add any number of specimen requirement groups to a collection event. However, the same specimen requirement may be a part of only one specimen requirement group.  
Select *Save study*.