

Record the collection date

Select the checkbox next to a specimen you have collected, or you can let the system do this for you after you have entered the date. Select the Calendar icon to record date and time of collection.

Collected	Study	Specimen expected	Quantity	Collection	Status	Actions
<input checked="" type="checkbox"/>	10603	3ml Whole Blood (Green Top)	3.0 ml	<input type="text"/>	Pending	
<input type="checkbox"/>	10603	5 ml Bone Marrow Aspirate (Green Top)	5.0 ml	<input type="text"/>	Pending	

Select the checkbox next to the specimen you have collected and system will populate the collection date with current day's date and time. You may, however, change the date and time later. Alternatively, select the calendar icon beside the collection date field after which system displays a calendar where you may select the date and time of collection.

SPECIMEN CHECKLIST FOR PATIENT 5000 (A,A,A) FOR STUDY 10603

Save Add to shipments

Pre-Registration Remission Induction Therapy Remission Consolidation Therapy Continuation Therapy Remission Relapse

Collected	Study	Specimen expected	Quantity	Collection	Status	Actions
<input type="checkbox"/>	10603	3ml Whole Blood (Green Top)	3.0 ml	<input type="text"/>	Pending	
<input type="checkbox"/>	10603	5 ml Bone Marrow Aspirate (Green Top)	5.0 ml	<input type="text"/>	Pending	

Now Done

Save Add to shipments

Record *Collection date* by selecting the desired date from calendar. To select time of collection, drag the sliders to adjust hours and minutes in time. If you have collected the specimen right now, select *Now* in the calendar and the *Done*.

Following is a screenshot indicating how the checklist view looks once *Collection date* is recorded.

Collected	Study	Specimen expected	Quantity	Collection	Status	Actions
<input checked="" type="checkbox"/>	10603	3ml Whole Blood (Green Top)	3.0 ml	06/07/2012 16:34	Pending	
<input type="checkbox"/>	10603	5 ml Bone Marrow Aspirate (Green Top)	5.0 ml	<input type="text"/>	Pending	

Once date and time of collection are entered, system selects the checkbox to indicate you have captured the collection date of that specimen. However, status of the specimen will still be *Pending* until You select *Save* located on top/bottom right corner to mark the specimen as collected. Status of the specimen is updated from *Pending* to *Collected* after you save Specimen checklist of that patient

Placebo : Cycle 1-Day 21									
<input checked="" type="checkbox"/>	10603	3ml Whole Blood (Green Top)	3.0 ml	06/07/2012 16:34	14	Collected			
<input type="checkbox"/>	10603	5 ml Bone Marrow Aspirate (Green Top)	5.0 ml		14	Pending			